

Power Drive Controls

1846 Santo Domingo
Camarillo, CA 93012
U. S. A.

www.pdcontrol.com

email: dashengpan@yahoo.com

Operation Notes for Red 650 nm SLM Laser

Module R650C

(Rev. 5-24-23)

1. Refer to the included **Panel Layouts** drawing for the module's operation functions
2. Plug the included **+5VDC wall power supply** into the **power jack**.
3. Turn the **power switch ON**.
4. The **Laser Enable LED** will turn **Red** then switch to **Green** after a few seconds.
5. Once the “**Temp Good**” LED is stabilized on **Green**, press the “**Laser Enable**” pushbutton switch to enable the laser
6. The module **pre-tuned to SLM**, no user adjustments are required.

Warning: Avoid **direct exposure** to the laser beam.